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# **OLYMPUS Cleanliness Inspector CIX ASW**

Quick Setup Guide

## TECHNICAL SUPPORT

If you experience technical problems, please contact our Technical Support by e-mail or telephone (Monday to Friday, 9:00 am to 5:00 pm Central European Time).

HOTLINE: +49 251 798 00 6444

E-MAIL: support@olympus-sis.com

When contacting our Technical Support, please have the following information to hand:

- Software version
- System name
- License number
- Microsoft Windows version
- Operating system language

Use the [System Information] > [About] > [More System Information] command to find all of the information that you need.

## NOTE

The OLYMPUS Cleanliness Inspector Software (CIX-ASW) has already been installed and activated on your OLYMPUS Cleanliness Inspector System. The following information only applies when the software is being reinstalled.


This software uses a license protection scheme which requires Internet-based activation. We recommend that you connect your computer to the Internet before installing the OLYMPUS Cleanliness Inspector Software.

A version of Acrobat Reader has been installed on your system to enable you to view reports.

A version of Acrobat Reader is also available on the installation DVD via the [Install example data or tools] option.

## INSTALLATION


To install your software, you need to have a valid license certificate. This license certificate is provided to you in a separate envelope. Open the envelope and look for the license key (15 characters in the blue area).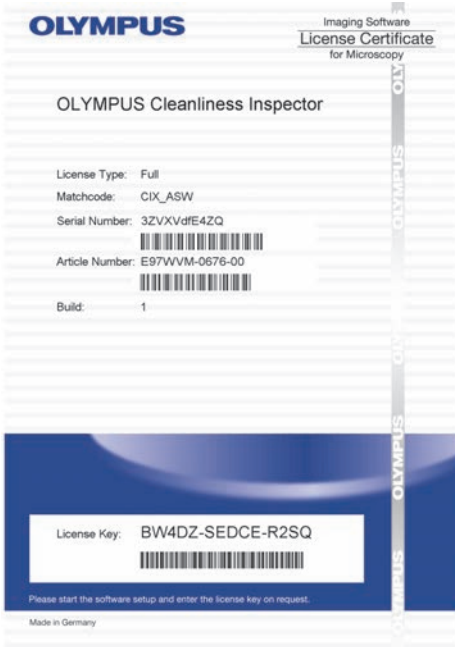

Insert the DL-DVD into your DL-DVD drive. The start dialog box of the installation program should be automatically displayed on your screen.

**Note:** It is possible that the autorun function of your operating system is deactivated. In this case, double click the “setup.exe” file located in the root directory of your DL-DVD drive.

In the first dialog box, select the [Install or maintain imaging software] option and click [Next >].

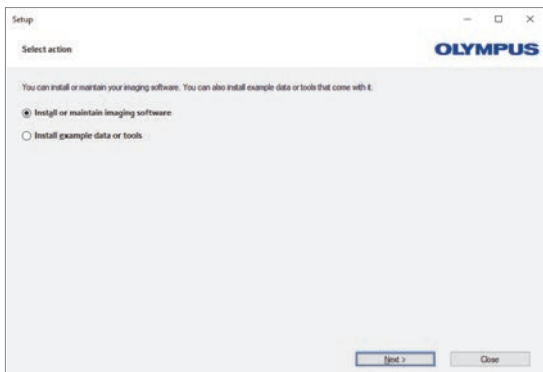The installation program's next dialog box asks you to enter your license key. Please enter all 15 characters into the required fields.

Now you can install and check your software straight away without having to activate it. Activation is compulsory and must be done within 30 days after the first installation.


## ACTIVATION

The license has to be activated within 30 days. Activation is not necessary for the 30 days demo license.


Activation requires a working Internet connection. The activation takes place in the background.

Click the [Contact Server] button to complete the activation of the OLYMPUS Cleanliness Inspector Software.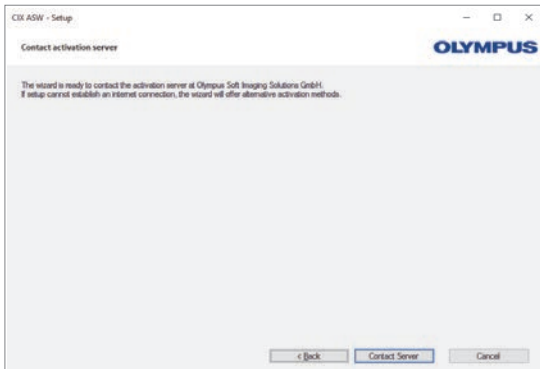


**Note:** If you are not connected to the Internet, but are able to exchange files with another computer that does have an Internet connection, please follow the steps in the next dialog box.

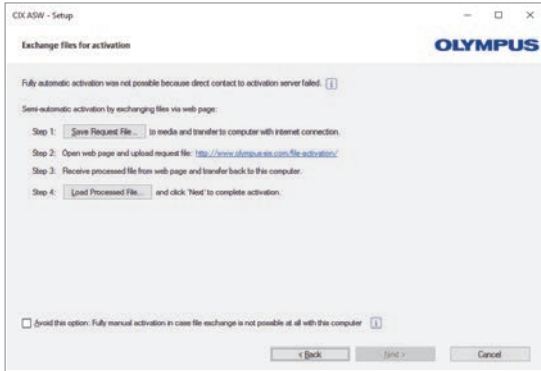The following website is used for the activation.

**Note:** If you are not connected to the Internet and are not able to exchange files with another computer that does have an Internet connection, you can activate the OLYMPUS Cleanliness Inspector Software by telephone. Call our Customer Service Center and follow the instructions below.

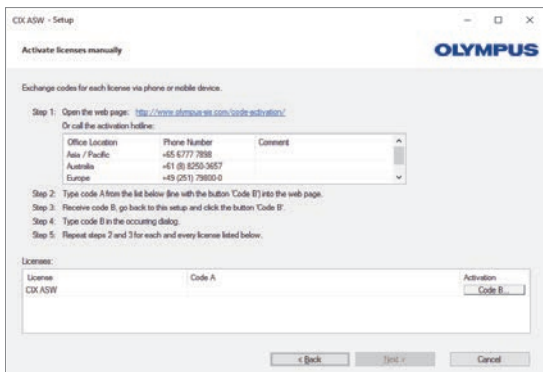The activation is now complete.

You can use your software on the computer for which it was activated, even if it doesn't have an Internet connection.

## DEACTIVATION

You must deactivate the OLYMPUS Cleanliness Inspector Software if you want to reinstall it on the same computer or install it on a different computer. Note the following advice. It forms part of the OLYMPUS Licensing conditions:

Deactivate your software license before you

- install and activate the same OLYMPUS Cleanliness Inspector Software on a different computer
- use a previously created restore point
- restore your system using a backup file
- replace hardware components on your computer
- restore or change your operating system.

Click the [System Information] > [Deactivate License] button on the software's start page. The installation program will lead you through the deactivation step-by-step. You can also start the deactivation process with the installation program on the setup DVD.

## SOFTWARE UPDATES

Please contact the local distributor from whom you acquired the product to inquire about software updates for the OLYMPUS Cleanliness Inspector Software (CIX ASW).

# **OLYMPUS Cleanliness Inspector CIX ASW**

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