


SWTLU-C

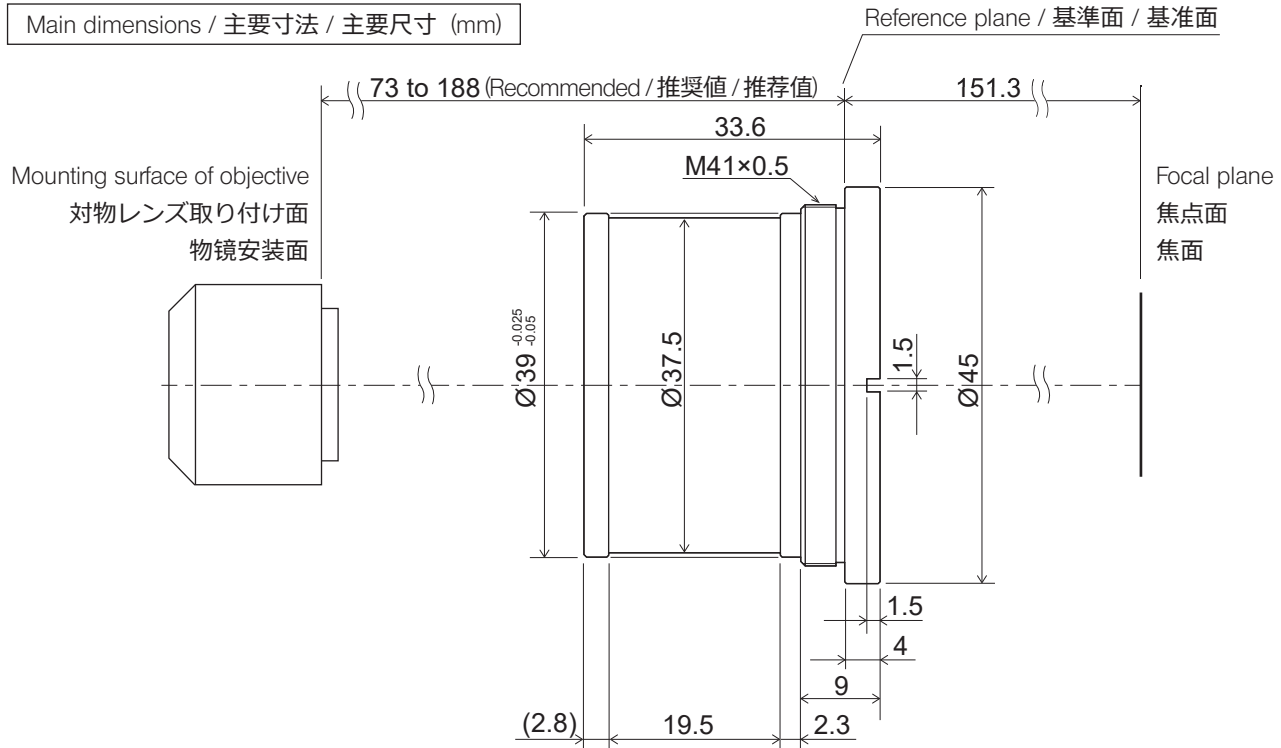

Precautions for use / 使用上の注意 / 使用注意事项

If the lens gets dirty, blow away dust with a commercially available blower and gently wipe the lens with a piece of cleaning paper (or clean gauze). Only when cleaning fingerprints and oil stains, slightly moisten a piece of cleaning paper with commercially available absolute alcohol and wipe the lens off with it.

The surface treatment for stray light prevention is applied to the lens frame of this product. Be sure not to wipe the lens frame when wiping the lens, otherwise the surface treatment film could be removed.


レンズが汚れた場合には、ほこりを市販のプロアーなどで吹き飛ばし、クリーニングペーパー（または洗いざらしの清潔なガーゼ）で軽く拭く程度にしてください。指紋や油脂類の汚れのみ、市販の無水アルコールをクリーニングペーパーにわずかに含ませて拭き取ってください。

この製品のレンズ枠には、迷光防止のための表面処理が施されています。レンズ枠を拭くと表面処理がはがれますので、レンズを拭くときにはレンズ枠を拭かないようにご注意ください。

如果在镜片上附着脏污时，请使用市售的气吹球等吹掉灰尘后，再用清洁纸（或洁净的纱布）轻轻擦拭镜片。

如果有指纹和油污时，请使用蘸上市售的无水酒精的清洁纸轻轻擦拭。

在本产品的镜片框上施加防止杂散光的表面处理。如果擦拭镜片框，表面处理将会剥落。擦拭镜片时切勿擦拭镜片框。

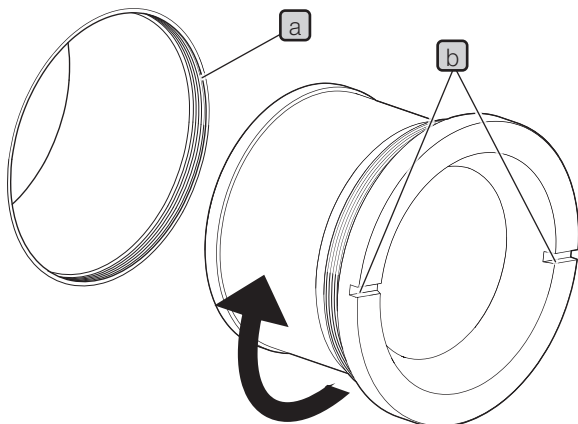

Installation / 取り付け / 安装

Screw the SWTLU-C into the M41X0.5 screw hole **a**. After the SWTLU-C is screwed in to the end, insert tools into the slots **b** (2 places) and rotate them clockwise to tighten the SWTLU-C securely.

SWTLU-CをM41X0.5のねじ穴**a**にねじ込みます。最後までねじ込んだら、すり割り**b**(2カ所)に工具を挿し込んで時計回りに回し、しっかりと固定します。

将SWTLU-C给M41X0.5的螺丝孔**a**拧进去。拧进到底后，将工具插入到产品狭槽**b**(两处)，正转工具以牢固固定SWTLU-C。

EVIDENT CORPORATION
株式会社エビデント

销售商名称 - 地址

仪景通光学科技(上海)有限公司

光学显微镜附件

总部：中国（上海）自由贸易试验区日耀北路199-9号102及302部位
北京分公司：北京市朝阳区酒仙桥路10号1幢一层102室
售后服务热线：400-969-0456