CODED 7 POSITION NOSEPIECE / CODED 6 POSITION NOSEPIECE

U-D7RES / U-D6RES



Connecting the coded revolving nosepiece (U-D7RES/U-D6RES) to the microscope frame (BX43 or BX53F2) makes it possible to use the light manager function of the microscope. Also, connecting the coded revolving nosepiece (U-D7RES/U-D6RES) to the microscope frame (BX63F) or the control box for coded function (U-CBS) makes it possible to read out the hole position information of the coded revolving nosepiece.

1 Cable connection

Connecting to the microscope frame

Lay out the cable by attaching the cable holders provided with the coded revolving nosepiece on the back side of the microscope frame as shown below.

- * If the cable holder positions are significantly deviated, the cable may not reach the connector. Attach the cable holders while making sure that the cable reaches the connector.

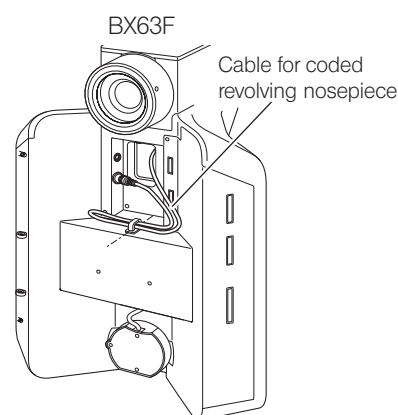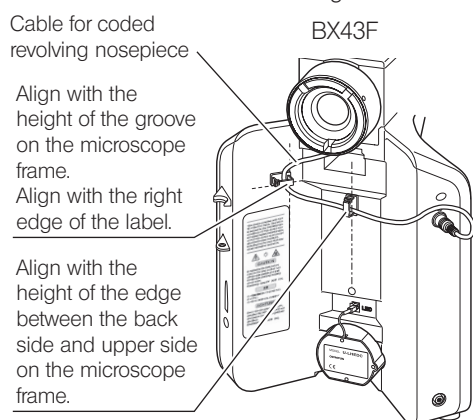If you connect to the microscope frame (BX53F2), refer to the instruction manual provided with BX53F2.

When connecting to the control box for coded function (U-CBS) or the interface for coded nosepiece (U-IFRES)

Refer to the instruction manual provided with U-CBS.

2 Operating procedures

For the operating procedures, refer to the instruction manuals provided with BX43F, BX53F2, BX63F and U-CBS.

- * When combining U-D6RES with BX63F, select [U-D7RES] in the revolving nosepiece settings on the touch panel controller provided with BX63F.

3 Specifications

| Item | Specifications | |
|---------|----------------|-------------------------------------|
| U-D7RES | Dimension | 119.5 (W) x 134.5 (D) x 47.2 (H) mm |
| | Weight | 1.1 kg |
| U-D6RES | Dimension | 107.5 (W) x 116 (D) x 48.2 (H) mm |
| | Weight | 0.7 kg |

Before disposing of this product, be sure to check the regulations and rules of your local government and follow them. Contact Olympus for any questions.

This product is a part of the system product which complies with the CE marking.
Please refer to the instruction manual of your system product for the safety instructions related to the CE marking.

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