# U-IFRES


# INSTRUCTIONS

Combining this product with the microscope frame (BX43F, BX53F2) and the control box for coded function (U-CBS) makes it possible to use the light manager function of the microscope frame and the function to readout the revolving nosepiece information to the digital camera/imaging software at the same time.

## 1 Cautions

- 1) For operating procedures of the microscope system attached with this product, refer to the instruction manuals provided with the microscope frame and U-CBS.
- 2) Before connecting the cables, make sure that the powers of the microscope frame and U-CBS are set to OFF.
- 3) Do not stretch or bend the cables with the excess force. Otherwise, they may be damaged. When connecting the cables, be sure to lay out the cables according to the instruction manual provided with U-CBS.
- 4) When carrying the microscope system, be careful so that the cables are not pinched under the microscope frame.
- 5) Before disposing of this product, be sure to check the regulations and rules of your local government and follow them. Contact Olympus for any questions.

## 2 Attaching the body of interface for coded nosepiece (U-IFRES)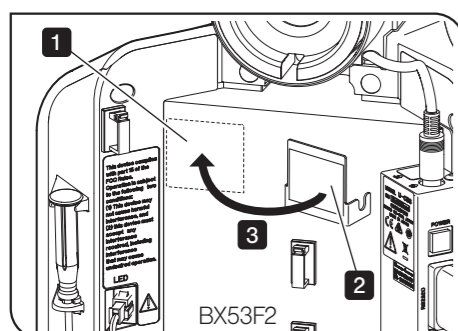


- 1 Clean the area to attach the fixing board on the back side of the microscope frame with a soft cloth.
- 2 Remove the protection sticker from the fixing board.
- 3 Attach the fixing board to the back side of the microscope frame firmly.

**TIP** If you remove the sticker of the fixing board attached once, the adhesive power may be weakened and U-IFRES may drop. Decide the position carefully before attaching to avoid the double attachment.

### When using BX53F2


Align the fixing board with the upper left of the back side of the microscope frame.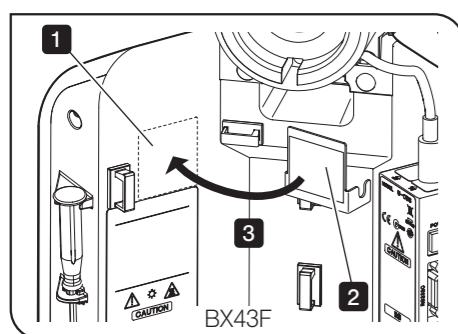

### When using BX43F

Align the fixing board with the upper right corner of the label on the back side of the microscope frame.


**TIP** Be careful not to overlap with the label.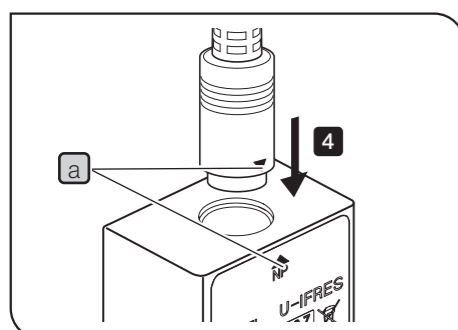

- 4 Connect the connector of the coded revolving nosepiece to U-IFRES.


**TIP**

- Insert the connector by matching ▼ marks **a**.
- For coded revolving nosepiece, refer to the instruction manual for U-D7RES/U-D6RES.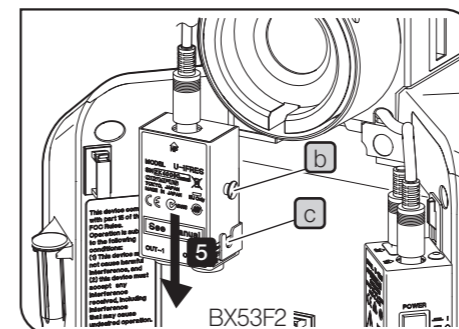

- 5 Attach U-IFRES by hooking the pin **b** of U-IFRES on the groove **c** of the fixing board.

**TIP** If you attach or remove the connector while U-IFRES is being hooked on the fixing board, the sticker of the fixing board may come off and U-IFRES may drop. When you attach or remove the connector, be sure to hold U-IFRES with your hands.

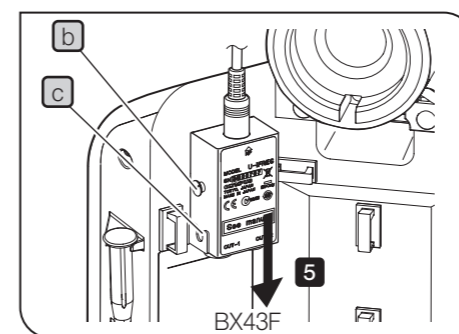## 3 Layout of the cable

Refer to the instruction manual provided with the U-CBS.

## 4 Specifications

Item	Specifications	
U-IFRES	Dimension	Body: 44 (W) x 25 (D) x 60 (H) mm Fixing board: 40 (W) x 19 (D) x 38 (H) mm
	Weight	0.15 kg
Operating environment	<ul style="list-style-type: none"> <li>• Indoor use</li> <li>• Altitude: Max. 2,000 meters</li> <li>• Ambient temperature: 5 to 40 °C (41 to 104 °F)</li> <li>• Maximum relative humidity: 80% for temperatures up to 31 °C (88 °F) (without condensation) In case of over 31 °C (88 °F), the relative humidity is decreased linearly through 70% at 34 °C (93 °F), 60% at 37 °C (99 °F), and to 50% at 40 °C (104 °F).</li> <li>• Supply voltage fluctuations: Not to exceed ±10% of the normal voltage.</li> <li>• Pollution degree 2 (in accordance with IEC60664-1)</li> <li>• Installation/Overtoltage category: II (in accordance with IEC60664-1)</li> </ul>	

### Repair request

If the abnormality or defect is found with the product, contact Us for assistance. At that time, please tell them the following information as well.

- Product name and abbreviation (Example: interface for coded nosepiece U-IFRES)
- Serial No. (SN)
- Problem

This product is a part of the system product which complies with the CE marking. Please refer to the instruction manual of your system product for the safety instructions related to the CE marking.

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