
Extendable eyepoint adjuster SZX2-EEPA INSTRUCTIONS

This product is the "extendable eyepoint adjuster" used in combination with the stereo microscope series SZX7, SZX10 or SZX16. For usage not described in this instruction manual, refer to the instruction manual for your microscope.

1 Handling precautions

- This product is a precision instrument. Handle it with care and avoid subjecting it to a sudden or severe impact.
- Never disassemble any part of the product. Otherwise, failure could be caused.
- For cleaning, storage and operating environment of this product, refer to the instruction manual for your microscope.
- Before disposing of this product, be sure to follow the regulations and rules of your local government.
- Due to performance limitations, this product cannot be used in combination with the following products:
 - Intermediate attachments: SZX-SDO2
 - Fluorescence units: SZX2-RFA16 and SZX2-RFA16A
 - Stands: SZ2-STU1 and SZ2-STU3
 - Camera: DP80
- Do not use this product with the following product combinations. Otherwise the microscope may be tilted or fall over.
 - LED transmitted light illumination base (SZX2-ILLTQ/ILLTS) + Optional column (SZH-P400/SZH-P600) + Fluorescence units

2 Adjusting eyepoint height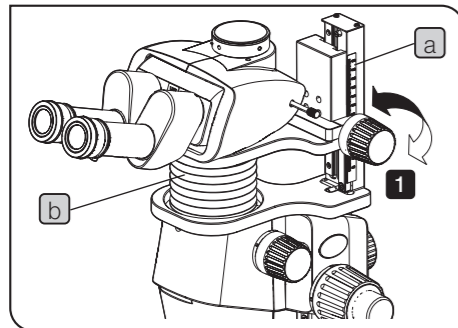

- 1 Turn the eyepoint height adjustment knob to move the movable part up and down.

The eyepoint is raised/lowered by turning the knob clockwise/counterclockwise.

TIP The eyepoint height (difference from the height without the eyepoint adjuster) can be checked by the scale **a** located on the right side. If you take note of your optimum eyepoint height, it is convenient to read just the height easily next time.

NOTE • The eyepoint height should be set within the range from 30 to 150 mm of the scale **a**. If you use it outside of the range (close to both ends of the height adjustment mechanism), the observation image may be moved remarkably. Also, if the movable part is pushed to the stoppers at both ends of the height adjustment mechanism with a strong force, the eyepoint height adjustment knob may rattle.


• Do not apply excessive force to the bellows **b** covering between the movable part and the fixed part of the eyepoint adjuster. Otherwise the bellows may be deformed.

3 Specifications

Item	Specifications
Eyepoint height adjustment range	30 to 150 mm (difference from the height without the eyepoint adjuster) Scale indicator is available
Total loaded weight limit	5 kg
Dimension	227(W) x 198(D) x 194(H) mm
Weight	2.1 kg


4 Attaching the eyepoint adjuster and the observation tube

Use the Allen screwdriver provided with the stereo microscope to attach the units.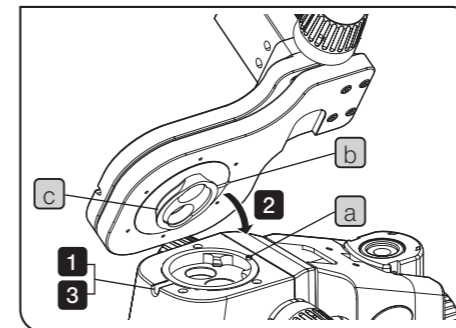

- 1 Loosen the clamping screw of the observation tube mount of the zoom microscope body.
- 2 Aligning the positioning pin **a** of the observation tube mount to the positioning groove **b** on the eyepoint adjuster, insert the male dovetail **c** at the bottom of the eyepoint adjuster into the observation tube mount.
- 3 While holding the eyepoint adjuster firmly in one hand, tighten the clamping screw of the observation tube mount securely.

NOTE If the eyepoint adjuster is released before it is clamped securely, it may be removed from the zoom microscope body and fall down.

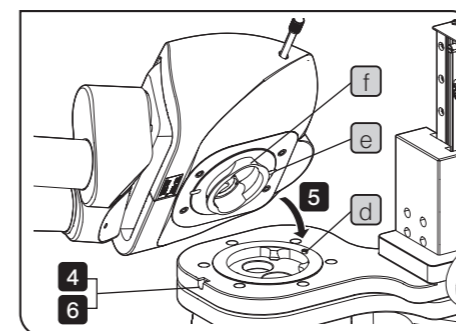- 4 Loosen the clamping screw of the observation tube mount of the eyepoint adjuster.
- 5 Aligning the positioning pin **d** of the observation tube mount to the positioning groove **e** on the observation tube, insert the male dovetail **f** at the bottom of the observation tube into the observation tube mount.
- 6 While holding the observation tube firmly in one hand, tighten the clamping screw of the observation tube mount securely.

NOTE If the observation tube is released before it is clamped securely, it may be removed from the eyepoint adjuster and fall down.

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