

The U-LHAD is a set of adapters for use in attaching the U-DP dual port onto the lamp housing in order to combine it with the mercury burner and other light source.

1 APPEARANCE, NOMENCLATURE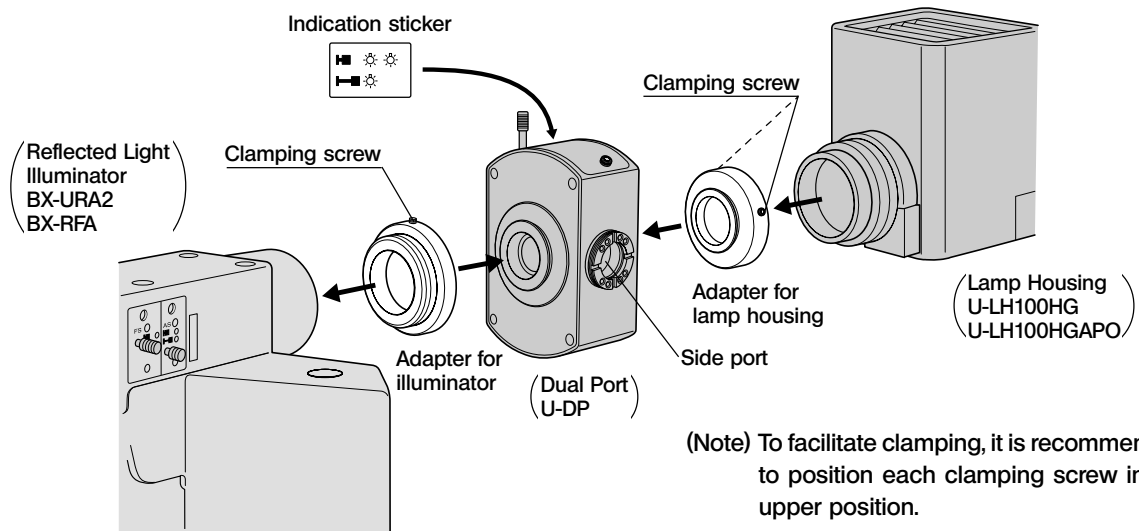

2 ASSEMBLY

Dimensions of Side Port

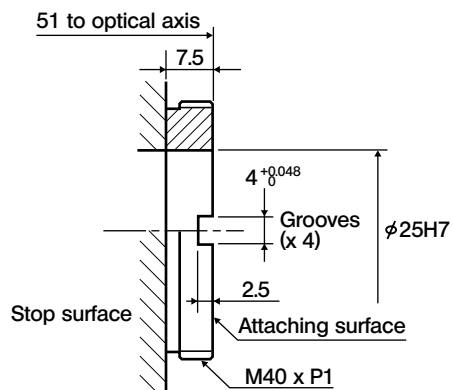⊙ Connect modules as shown in the above illustration.

Attach the provided indication stickers above the indications on the U-DP.

⊙ The adapter for use in attaching a lamp housing or similar module onto the side port can be fabricated on custom order.

If you want to fabricate such an adapter by yourself, refer to the dimensions shown in the following illustration. When a parallel beam is incident to the port, approximately the entire specimen will be illuminated. For the detailed optical data, please consult your dealer.

3 RESTRICTIONS

- When the U-LHAD is used, the field number of the illumination system becomes 11.
- The weight of the light source, etc. attached on the side port should be no more than 1 kg. (If the weight is 1 kg or more, attach an additional support.)
- The U-DP1XC can be attached to the side port but the observation image cannot be obtained.

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