OLYMPUS

Optical Microscope Accessory

Safety and Warning Information Digital Microscope Cameras

This flyer applies to the following cameras: DP23, DP23M, DP28, LC35, SC180. You can download the complete instruction manual of your camera and available software from the [Support] section of the following websites. Choose [Software Downloads] or [Instruction Manual Download].

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Digital Microscope Camera

Instruction manual

Before using the camera, download the complete instruction manual. In order to be able to read the instruction manual, software that can read PDF files must be installed on your computer. Make sure that you have access to the complete instruction manual when you are using the camera.

Safety

The camera has been developed and produced in conformity with recognized safety regulations. However, it is not possible to entirely rule out possible hazards for the user, damage to the camera or microscope, or impairments of the camera's functionality. For this reason, always read and follow the general safety instructions and warnings before and during installation.

⚠ WARNING - Avoiding personal injury

Electric shock due to damaged cable

When a cable is frayed or damaged, a person touching it risks receiving an electric shock. Shut off the hardware and replace the cable immediately.

Poisoning due to toxic gases

In the event of a fire, the material of the camera can give off toxic gases. When these are inhaled, it can cause irritation or damage to the respiratory tract. Wear

respiratory protection when fighting the fire.

⚠ CAUTION - Avoiding personal injury

Tripping hazard

Cables laid across a room can be a dangerous tripping hazard. This can cause injuries to people and damage to equipment. Wherever possible, lay cables along the wall or behind furniture. Fix and label cables that run across the room.

ATTENTION - Avoiding device damage

Never open the camera

The camera's CMOS chip is extremely sensitive. Electrostatic discharge can permanently damage the camera. Therefore, never open the camera housing.

Use only the supplied USB camera cable

Operate the camera only with the supplied USB camera cable. Only then will the functionality of the camera and compliance with the EMC basic standards be ensured.

Don't touch the protective glass

The protective glass over the CMOS chip is extremely sensitive and must not come into contact with bare hands or any other objects, otherwise the acquired images will be impaired by fingerprints or scratches.

Digital Microscope Camera

Avoid vibration shocks

Mechanical shocks and strong vibrations can damage the camera. Don't drop the camera during assembly. Make sure the camera is securely mounted on the microscope. When the camera is not in use, store it in a safe place.

Comply with operating and storage conditions

Exposure to high temperatures or humidity can damage component parts of the camera. Always observe the operating and storage conditions.

Explanation of the symbols used

WARNING

Indicates a potentially hazardous situation which, if not avoided, may result in major injury.

CAUTION

Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.

ATTENTION

Indicates a situation which, if not avoided, may damage the equipment or other property.

Contacting Customer Service

Please contact your local Olympus support if you have any questions about the product. More information can be found on this website:

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EVIDENT Technology Center Europe GmbH
Evident Europe GmbH
Evident Scientific, Inc.
48 Woerd Ave Waltham, MA 02453, U.S.A.
Evident Korea Co. Ltd.
Evident Scientific Singapore Pte. Ltd.