## How to use U-D5BDREMC-VA3

- 1 Slowly remove the sticker on the hole you want to use with your precision screwdriver, etc.
- 2 Insert the fitting ( $\phi 4$  or  $\phi 6$ ) provided with the motorized nosepiece into the hole you want to use and tighten it with your Allen wrench.

- TIP**
- Do not overtighten the fitting. Overtightening the fitting will damage it.
  - Do not remove screws, etc. near the hole you want to attach the fitting.

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**NOTE:** This product has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the product is operated in a commercial environment. This product generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this product in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

**FCC WARNING:** Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the product.

**FCC Supplier's Declaration of Conformity**

Hereby declares that the product

Product name: Optical Microscope Accessory

Model Number: U-D5BDREMC-VA

Conforms to the following specifications:

FCC Part 15, Subpart B, Section 15.107 and Section 15.109

Supplementary Information:

This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

Responsible Party Name: Evident Scientific, Inc.

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This product is applied with the requirements of standard IEC/EN61326-1 concerning electromagnetic compatibility.

- Emission: Class A
- Immunity: Industrial electromagnetic environment

Emissions exceeding the level required by aforementioned standards may occur if this product is electrically connected to other equipment.

Some interference may occur if this product is used in residential environments.

In accordance with European Directive on Waste Electrical and Electronic Equipment, this symbol indicates that the product must not be disposed of as unsorted municipal waste, but should be collected separately.

Refer to your local Evident distributor in EU for return and/or collection systems available in your country.

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Address: 48 Woerd Ave Waltham, MA 02453, U.S.A.