## About Attachment of the zoom head

For the installation dimensions when incorporating the zoom head into your system, refer to the following dimensions for the mounting stand; provide a boss hole ( $\Phi 20$ ), M6 Tap (3 locations), and the butting pin. Make sure that there is no space between the zoom head and the stand. Otherwise, dirt/dust or water may enter, and the product may be damaged.



The cover is provided on the back of the zoom head in order to prevent the static electricity. Do not remove the cover. Be careful not to touch the board.

Use the zoom head tilt within the range of  $\pm 90^\circ$ . Rotate the revolver in the upright position.


If the product is installed in your system, the manufacturer provides no guarantee of calibration or measurement accuracy. Make sure that the product with zoom head combined complies with regulations required by the location (country) where you use this product.

Considering the vertical movement range of the zoom head, leave a space of at least 285 mm above the center of the boss hole. (The zoom head will rise 53mm from the position where it was attached at initialization.)

(Unit: mm)  
The tolerance is a recommended value.

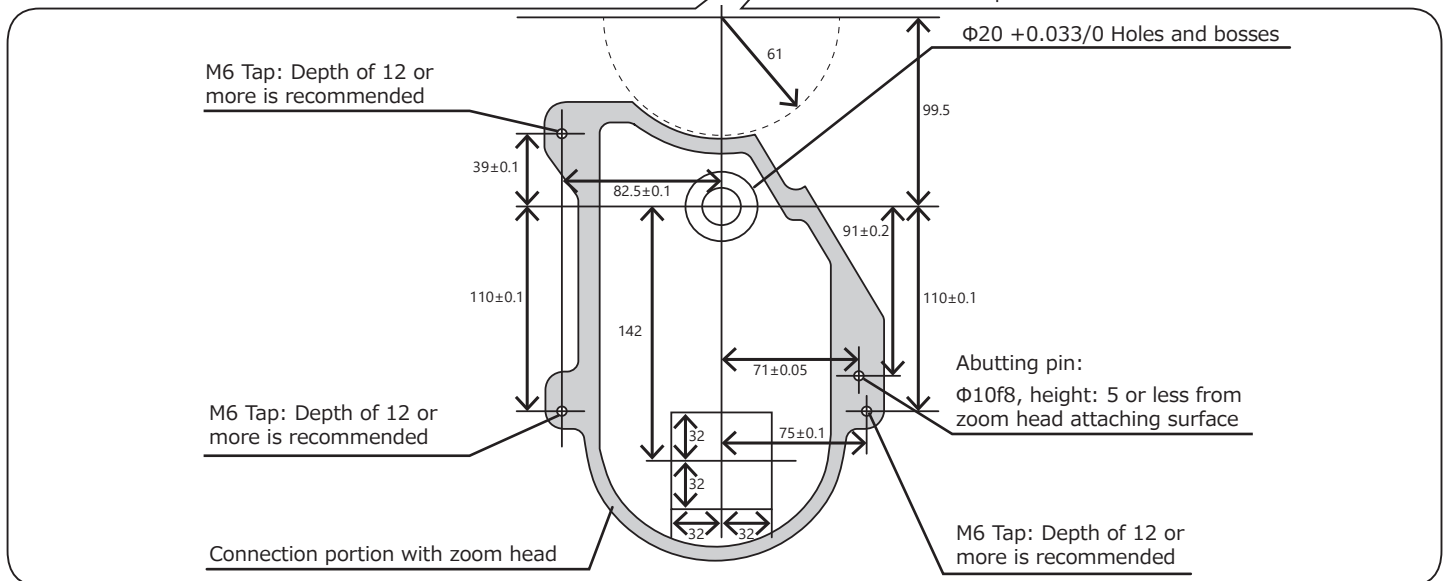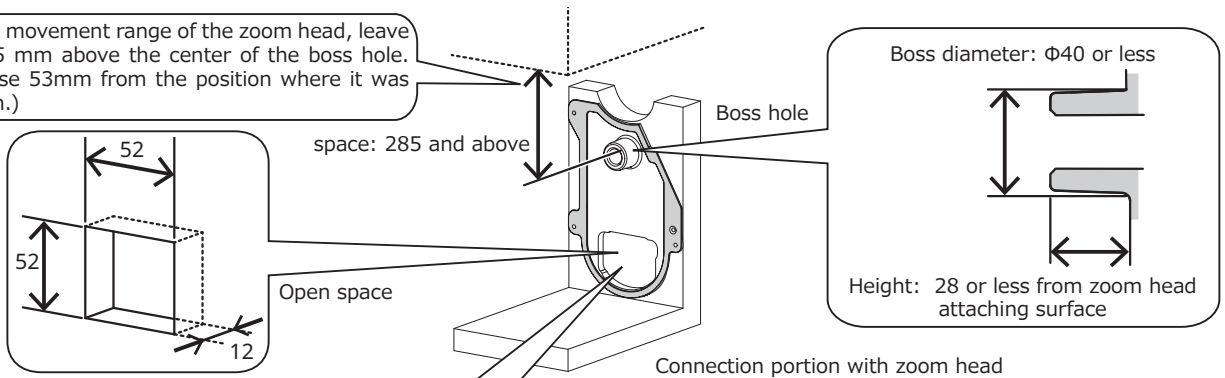### Zoom head Range of motion
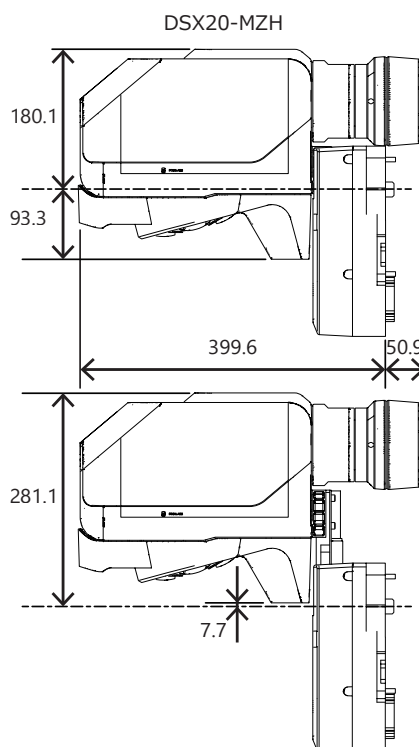
## Evident Corporation

6666 Inatomi, Tatsuno-machi, Kamiina-gun, Nagano 399-0495, Japan

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Evident Customer Information Center

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[https://www.olympus-lifescience.com/  
support/service/](https://www.olympus-lifescience.com/support/service/)

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[https://www.olympus-ims.com/  
service-and-support/service-centers/](https://www.olympus-ims.com/service-and-support/service-centers/)

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<https://www.olympus-ims.com>